

Date: 12 Jan 2026

Revision: 1.0

Lysis plate pickup and Xpress-seq plate submission guidelines

Please read and follow these guidelines carefully to ensure optimal results and smooth handling of your Xpress-seq samples.

1. Pick up lysis plates.

Please visit our webpage (www.xpress-genomics.com/booking/) to book a time slot for picking up/shipping lysis plates. The shelf life of our lysis plates are 12 months when stored in -20°C.

2. Invoice.

Once you have received the lysis plates we will invoice you 500 SEK / 45 Euro per plate. The pre-paid cost of lysis plates will be deducted from the final invoice of the Xpress-seq plate sequencing service.

3. FACS-sort.

Fill as many lysis plates as you wish with your sample/s.

4. Submit Xpress-seq plates.

Please visit our webpage (www.xpress-genomics.com/booking/) to book a time slot for handing in/shipping the FACS-sorted plates. Alternatively, you can ship plates to us using courier (eg. DHL Express, FedEx, UPS). Always make sure to transfer the plates fully buried in dry ice.

Important: We only accept plate submissions accompanied by correctly filled submission forms.

Xpress Genomics AB

R8:00, Cancercentrum Karolinska
Visionsgatan 56
171 64 Solna
Sweden

Delivery Contact Phone: +46 760079794

Delivery Contact Email: lab@xpress-genomics.com